



Excitation Filter Assembly P/N 10-150:

Unscrew the nuts for the plate and remove it. Sandwich the filter between the plates and replace the nuts. Care should be taken to not get finger prints / dust on the filter. If the filter has a mirror make sure the mirrored side is the side that shows in the picture. See next slide for install.



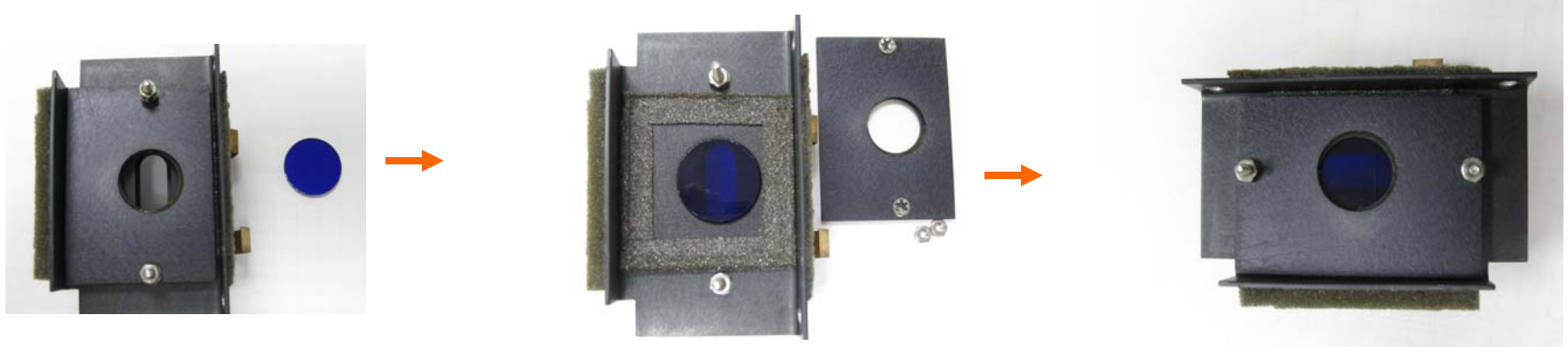


Emission Filter Assembly P/N 1000-260 and Install:


Set the filter in the holder and secure in place with o-ring or clip provided. Mirror side up. Remove front cover, lamp and old filters. Place the filter holding in the orientation shown. *It may work best to lay the 10-AU down on it's back for better access. Make sure to remove fingerprints / dust.

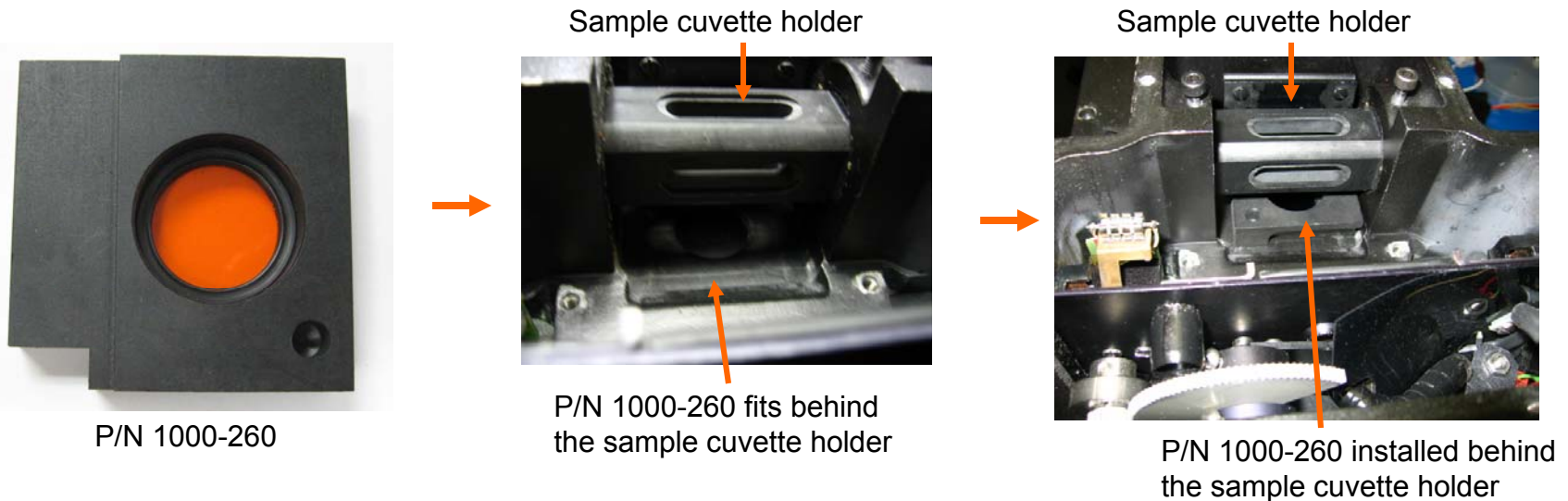

Install the Excitation Filter Assembly as shown. Secure with screws and replace lamp.

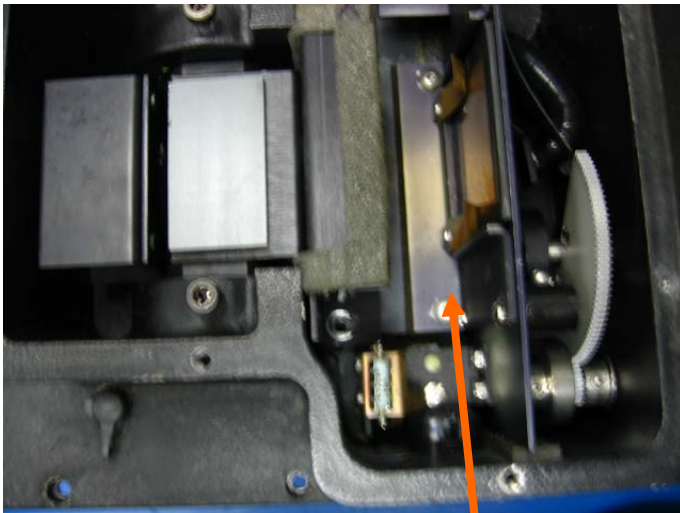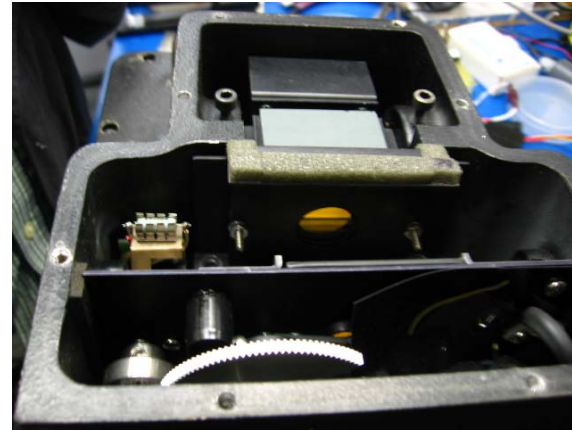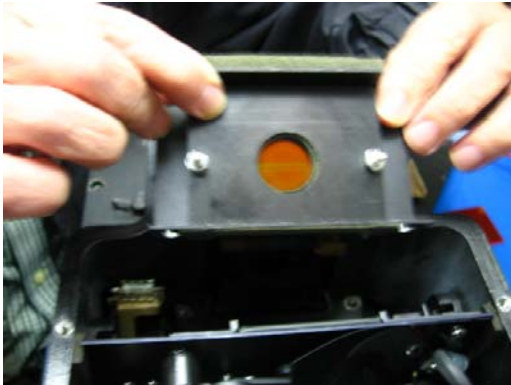

Lamp goes here.

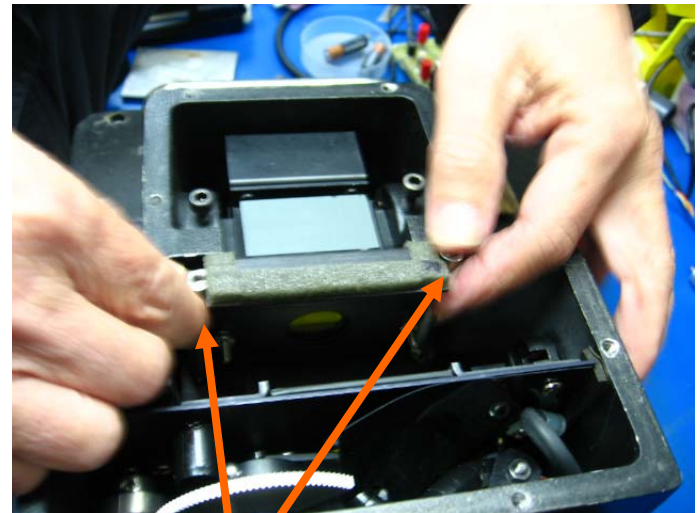

Long screws on each side of holder.