


CAUTION: MAKE SURE THAT THE DATABANK BATTERY IS FULLY CHARGED BEFORE ATTEMPTING UPGRADE.


To install or upgrade your DataBank instrument with the latest firmware version 3.15 and software version 3.15, you must first uninstall your current version of the DataBank GUI Software using the 'Add/Remove Programs' function found in your computer's control panel.

After removing the DataBank GUI Software from the programs list follow steps 1-11

1. Download the "DataBank Software Ver. 3.15" zipped folder from the following link:
<http://docs.turnerdesigns.com/Software/DataBank-Software-Ver.-3.15.zip>
2. Unzip and open the folder, then double-click the 'setup.exe' icon to re-install DataBank GUI Software Version 3.15.
3. After the installation has completed, connect your DataBank to your computer using the DataBank interface cable
4. Turn on your DataBank instrument
5. Begin the DataBank GUI Software program which can be found in the C:\Program Files\TurnerGUI folder. At this point the DataBank GUI Software Version 3.15 will give you an upgrade notice

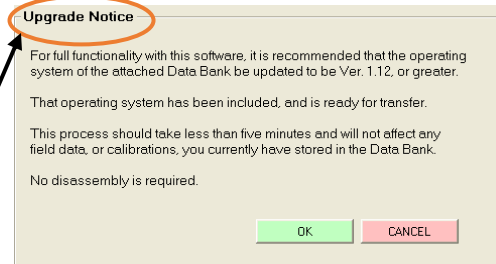

6. Click 'OK' to allow the upgrade to continue
7. Click 'Browse' and navigate to C:\Program Files\Turner GUI
8. Select the file titled 'DB315.HEX' and the following window will appear

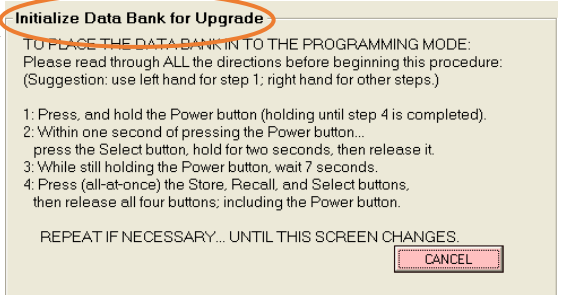

9. Follow the instructions in the window to enter programming mode
10. When initialization is triggered, a new window will appear indicating programming mode has been entered and the DataBank will begin uploading the new version of firmware
11. When firmware upload has completed, you will be instructed to reset the DataBank (follow onscreen instructions)

When DataBank is powered after the reset, the new software version 3.15 and firmware version 3.15 will be displayed in the main DataBank GUI Software window as shown below.

Software Version 3.15

Firmware Version 3.15

