

## 10-AU Calibration for Chlorophyll



### **3 Main Steps**

1. Setting the basic operating level
2. Blanking the instrument
3. Calibration with a liquid primary standard that has a known concentration

## #1 Setting the Basic Operating Level

This procedure should be performed the 1<sup>st</sup> time you use the instrument, or any time you switch to a different application, change filters or lamps.

1. Turn on the 10AU and allow the instrument to warm up for at least 30 minutes.



**2. Table 1 is an example of how set the sensitivity of the 10-AU. Users should determine the appropriate range and %FS for their calibration standard.**

Chlorophyll (ug/l)	Rhodamine WT (ppb)	Range	% FS (+/-5)
240 - 180	100 - 75	High	80%
180 - 160	75 - 65	High	70%
160 - 130	65 - 55	High	60%
130 - 110	55 - 45	High	50%
110 - 80	45 - 35	High	40%
80 - 60	35 - 25	High	30%
60 - 40	25 - 15	High	20%
24 - 18	10 - 7.5	Medium	80%
18 - 16	7.5 - 6.5	Medium	70%
16 - 13	6.5 - 5.5	Medium	60%
13 - 11	5.5 - 4.5	Medium	50%
11 - 8.0	4.5 - 3.5	Medium	40%
8.0 - 6.0	3.5 - 2.5	Medium	30%
6.0 - 4.0	2.5 - 1.5	Medium	20%
2.4 - 1.8	1.0 - 0.75	Low	80%
1.8 - 1.6	0.75 - 0.65	Low	70%
1.6 - 1.3	0.65 - 0.55	Low	60%
1.3 - 1.1	0.55 - 0.45	Low	50%
1.1 - 0.8	0.45 - 0.35	Low	40%
0.8 - 0.6	0.35 - 0.25	Low	30%
0.6 - 0.4	0.25 - 0.15	Low	20%


Table 1 is an example of how to set the scale for a given linear range of the 10-AU.

For this example we are using a **20µg/L** standard.

**3. Prepare your standards and acetone blank in test tubes so you are ready to set the sensitivity and calibrate the instrument.**


**Be sure to clear past calibrations on screen # 2.6**

4. Access screen 2.43 and set range control to manual and range to the appropriate level for your standard. For our example the range should be set to MED on Screen 2.42.

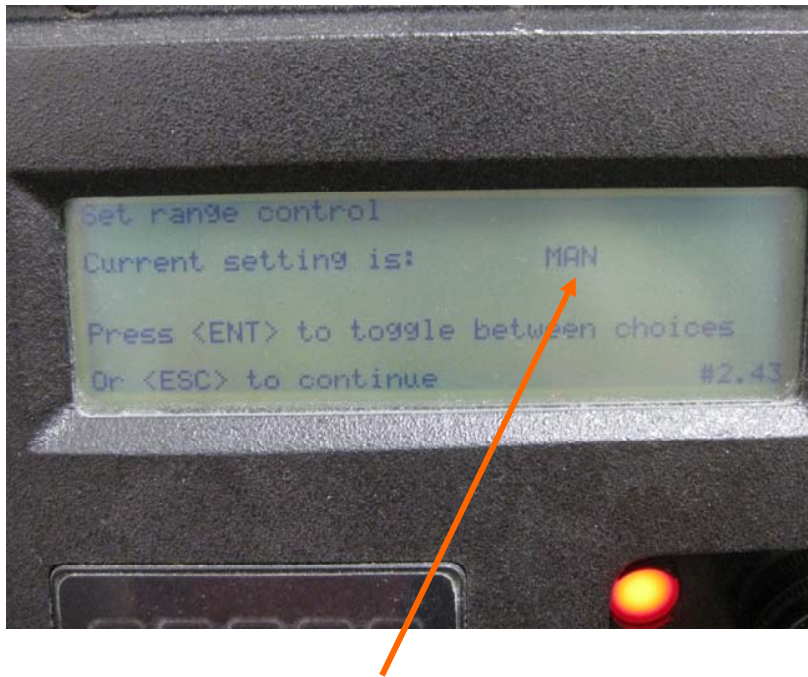

Change to MAN

#### 4. Access Screen 3.2

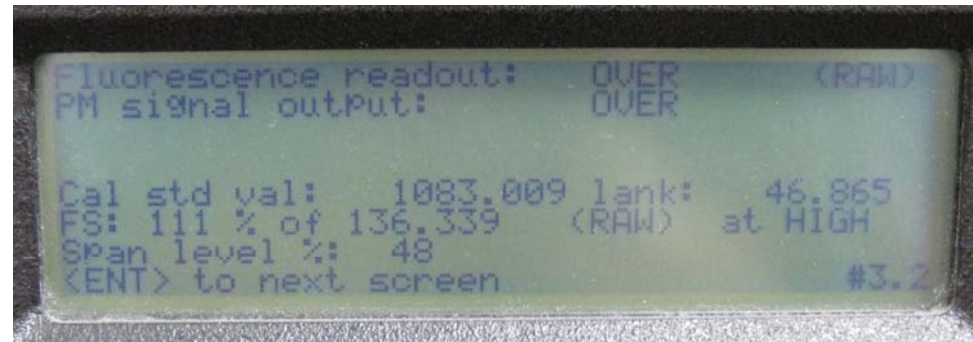

#### 5. Loosen the sensitivity locking screw using allen wrench.








8. When the %FS reaches the designated value, tighten the sensitivity locking screw. For our example this is  $\approx 80\%$  of medium



The Basic Operating Level is now set, and you can remove the standard for the 10-AU.

Setting the Basic Operating Level (aka Sensitivity) of your 10-AU is now complete.

You can access Screen 2.43 again and return range control to AUTO and begin calibration.

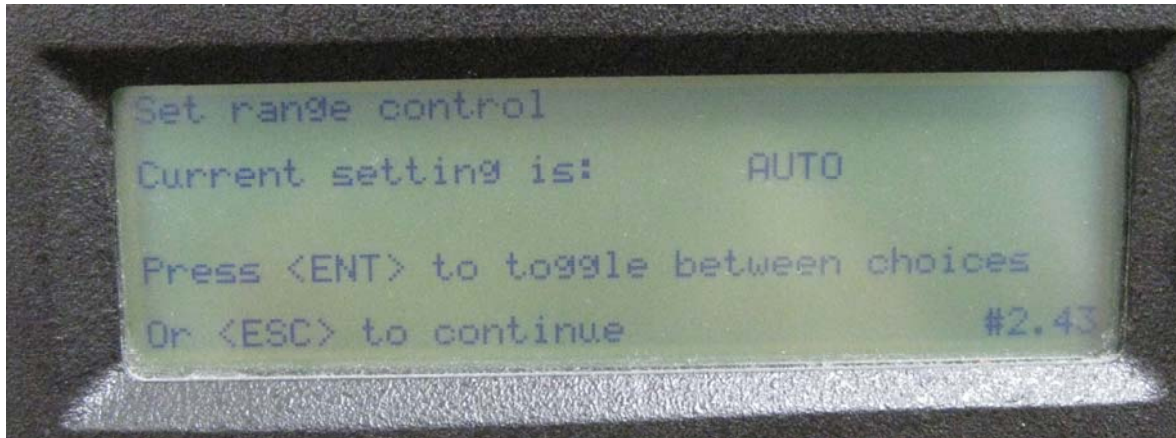

## #2 Blanking the Instrument

Now that the sensitivity is set, you should blank the instrument.

### 1. Access screen 2.1



### 2. Insert and cover your blank (acetone). Press 1 to Run blank.



**\*You can also choose to have the 10-AU subtract the blank signal from the sample readings on screen 2.12.**

Wait for the readings to stabilize and press “0”.

Wait for 15 seconds for the instrument to blank.


If Blanking is successful you will receive a  
“Blank Finished” message



## #3 Calibration with Your Standard

Now that the blank is set, you should calibrate with your standard.

### 1. Access screen 2.0

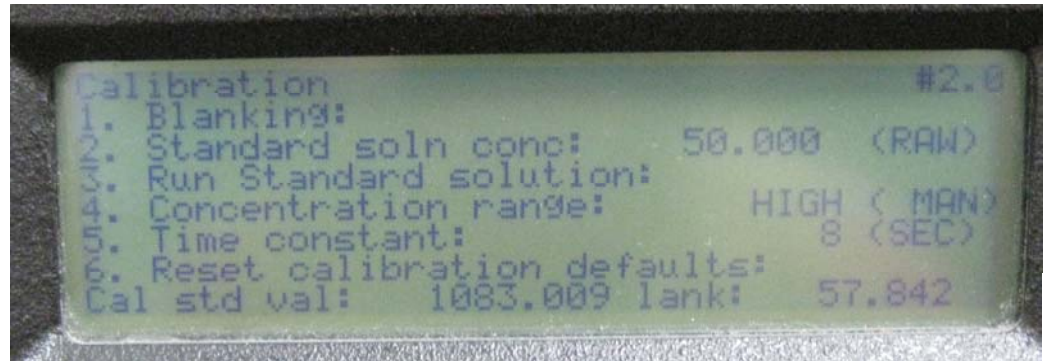


### 2. Press <2> to access screen 2.2 and input the actual concentration of your standard. Return to screen 2.0

### 3. Insert and cover your standard.



4. Press 3 to run standard. Wait for the reading to stabilize.

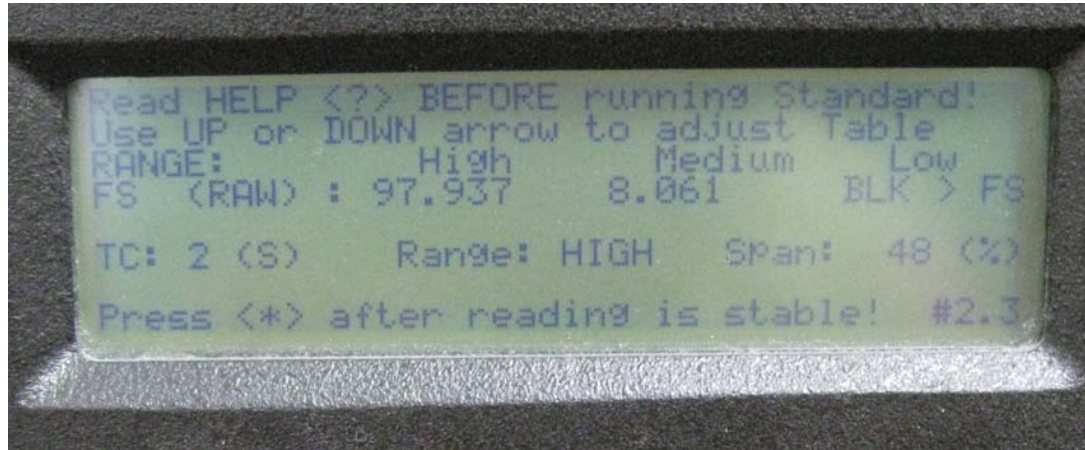

Press the <\*> key and the calibration point is set.  
Wait 15 seconds and you should receive a “FINISHED”  
message.




You can test/check your calibration by running your standards and other samples or standards.

I ran the 50 $\mu$ g/L and got:



I ran the blank and got:

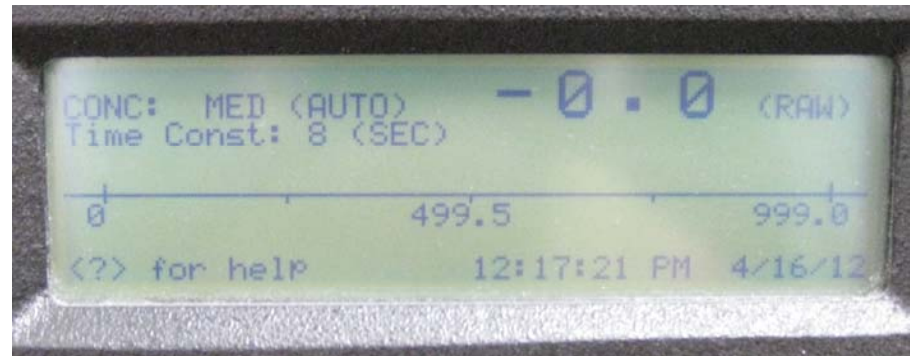

I ran a 10 $\mu$ g/L standard and got:

